



When product integrity counts...  
Manufacturers of laboratory calibrated temperature and humidity data loggers, real-time monitoring and cold chain solutions



## Instructions for installation of TRW 6.3 software

- Admin rights are required to install software
  - Do not plug in reader until software is installed.
1. Go to Temprecord website: <https://temprecord.com/our-software/temprecord-windows-software/>
  2. Copy software from Downloads folder to PC
  3. Right click trw-set up exe and “extract all”
  4. Double click on set up file or EXE file and follow installation prompts.
  5. Plug in reader
  6. Re-booting maybe required
  7. Open software
  8. Insert logger into reader (with logger facing up and hole exposed)
  9. Click space bar to access logger (The COM-PORT may need to be chosen)
  10. **Go to Parameters to enable use of the button on the logger**



For loggers that look like these, DO NOT use TRW software. For Mon-T2 data loggers please use TempWrite software at: <https://temprecord.com/our-software/tempwrite/>